

## Mouse anti-Epithelial Specific Antigen (Ep-CAM)

Cat. No.: BMS048 (16 ml Ready-to-use)

### Instructions for use

#### Intended Use

This antibody is designed for the specific localization of Epithelial Specific Antigen (Ep-CAM) in formalin-fixed, paraffin-embedded tissue sections.

Anti-Ep-CAM antibody is intended for research use only.

#### Specifications

<b>Specificity:</b>	Epithelial Specific Antigen (Ep-CAM)
<b>Immunogen:</b>	MCF-7 cells (human breast carcinoma cell line)
<b>Clone:</b>	Ber-EP4
<b>Isotype:</b>	Mouse IgG1 kappa
<b>Species reactivity:</b>	Human +, others not tested

#### Summary and explanation

Ep-CAM/Epithelial Specific Antigen, also known as epithelial antigen and epithelial glycoprotein, is expressed in the cytoplasmic membrane and the cytoplasm of nearly all epithelial cells. However, there are some exceptions like most squamous epithelia, hepatocytes, renal proximal tubular cells, gastric parietal cells, and myoepithelial cells. Sometimes focal positivity is observed in mesothelial cells undergoing reactive changes.

Ep-Cam is found in the majority of adenocarcinomas (50 – 100 % according to diverse studies) and in neuroendocrine tumours including small cell carcinomas.

30% of all renal and liver cell carcinomas are positive for Ep-Cam. Also basal cell and basosquamous carcinoma are anti-Ber-EP4 positive in almost all cases whereas the majority of malignant mesotheliomas is negative for Ep-Cam. In an appropriate panel this fact can be used to discriminate between this tumour type and adenocarcinomas.

#### Reagent provided

Mouse monoclonal antibody in TBS with carrier protein and preservative for stabilisation in the following format:

**Ready-to-use:** 16 ml (Cat. No. BMS048)

#### Dilution of primary antibody

None

#### Storage and handling

The antibody should be stored at 2-8°C without further dilution.

If necessary, dilutions of the antibody should be done with a suitable antibody dilution buffer (e.g. ZUC025 from Zytomed Systems). The diluted antibody should be stored at 2-8°C after use. Stability of this working solution depends on various parameters and has to be confirmed by appropriate controls.

The antibody provided is stable until the expiry date indicated on the label, if stored at 2-8°C. Do not use product after the expiry date. Positive and negative controls should be run simultaneously with all specimens. If unexpected staining is observed which cannot be explained by variations in laboratory procedures and a problem with the antibody is suspected, contact Zytomed Systems' technical support or your local distributor.

#### Precautions

Use through qualified personnel only.

Wear protective clothing to avoid contact of reagents and specimens with eye, skin and mucous membranes. If reagents or specimens come in contact with sensitive area, wash with large amounts of water.

Microbial contamination of the reagent must be avoided, since otherwise non-specific staining may occur.

ProClin300 is used for stabilisation. A Material safety data sheet (MSDS) is available upon request.

## Staining procedure

Refer to the following table for conditions specifically recommended for this antibody. Also refer to detection system data sheets for guidance on specific staining protocols or other requirements.

Parameters	Zytomed Systems recommendations
*Pre-treatment:	Enzymatic or Heat Induced Epitope Retrieval (for example in Citrate buffer pH 6, ZUC028)
*Control tissue	Colon carcinoma
*Working dilution	None
*Incubation time	30 - 60 minutes

## Quality control

The recommended positive control tissue for this antibody is a colon carcinoma. We recommend carrying out a positive and a negative control with every staining run. Please refer to the instructions of the detection system for guidance on general quality control procedures.

## Troubleshooting

If you observe unusual staining or other deviations from the expected results please read these instructions carefully, refer to the instructions of the detection system for relevant information or contact your local distributor.

## Expected results

The antibody stains positive in the cytoplasmic membrane and the cytoplasm of epithelial cells in formalin-fixed, paraffin-embedded tissue. The interpretation of the results is solely the responsibility of the user. Any experimental result should be confirmed by a medically established diagnostic procedure.

## Limitations of the Procedure

Immunohistochemistry is a complex technique involving both histological and immunological detection methods. Tissue processing and handling prior to immunostaining, for example variations in fixation and embedding or the inherent nature of the tissue can cause inconsistent results (Nadji and Morales, 1983). Endogenous peroxidase, alkaline phosphatase or biotin may cause non-specific staining depending on the detection system used. Tissues containing Hepatitis B Surface Antigen (HBsAg) may give false positive results with HRP (horse radish peroxidase) detection systems (Omata *et al*, 1980). Inadequate counterstaining and mounting can influence the interpretation of the results.

Zytomed Systems warrants that the product will meet all requirements described from its shipping date until the expiry date is reached, if the product is stored and utilised as recommended. No additional guarantees can be given. Under no circumstances shall Zytomed System be liable for any damages arising out of the use of the reagent provided.

## Performance characteristics

Zytomed Systems has conducted studies to evaluate the performance of the antibody utilising a standard detection system. The product has been found to be sensitive and specific to the antigen of interest with minimal or no cross-reactivity.

## Bibliography

Ordóñez NG. Adv Anat Pathol 13:16-25, 2006  
Ordóñez NG. Mod Pathol 19:417-428, 2006  
Carella R *et al*. Am J Surg Pathol 25:43-50, 2001  
Ordóñez NG. Am J Clin Pathol 109:85-89, 1998






Latza U *et al*. J Clin Pathol 43:213-219, 1990  
Nadji M and Morales AR Ann N.Y. Acad Sci 420:134-9, 1983  
Omata M *et al*. Am J Clin Pathol 73(5): 626-32, 1980

September 24, 2013

Rev: A0913

Doc: DBE\_BMS048

Explanations of the symbols on the product label:

<b>REF</b>	Bestellnummer Catalog Number Reference du catalogue		Verwendbar bis Use By Utiliser jusqu'à		Gebrauchsanweisung beachten Consult Instructions for use Consulter les instructions d'utilisation
<b>LOT</b>	Chargenbezeichnung Batch Code Code du lot		Lagerungstemperatur Temperature Limitation Limites de température	<b>RUO</b>	Nur für Forschungszwecke For Research Use Only Pour la recherche uniquement
<b>IVD</b>	In vitro Diagnostikum In Vitro Diagnostic Medical Device Dispositif médical de diagnostic in vitro		Achtung Warning Attention		Hersteller / Manufacturer / Fabricant Zytomed Systems GmbH • Anhaltinerstraße 16 14163 Berlin, Germany • Tel: (+49) 30-804 984 990 www.zytomed-systems.com